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HLA-B27 RQ Kit Manual

Autumn 2025, Version 2.1

For Real-Time PCR Detection of HLA-B27
For Research Use Only

 24 (Cat# HLA-B27RQ24)

 48 (Cat# HLA-B27RQ48)

 96 (Cat# HLA-B27RQ96)

 NG-WI-ASL-43-201

RUO

NovinGene ParsVira

No. 5, Iranshahr St, Tehran, Iran 1581633336.

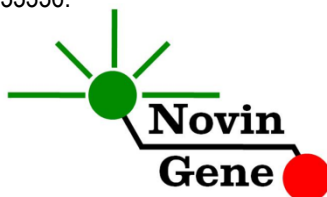
Table of Contents

1. Introduction	2
2. Intended Use	2
3. Background Information	2
4. Test Principle	3
5. Kit Contents.....	3
6. Packaging models	3
7. Storage and Stability	3
8. Product Use Limitations.....	3
9. Additionally Required Materials	4
10. General Precautions.....	4
11. Specimen, Storage and Transport.....	5
12. Interfering Substances	5
13. Internal Control.....	5
14. DNA Isolation	5
15. PCR Protocol	6
16. Devices and software	6
17. Programming Rotor-Gene	6
18. Programming StepOne.....	7
19. Programming Other Machines.....	8

20. Data Analysis: Rotor-Gene	8
21. Data Analysis: StepOne	11
22. Disposal Method.....	13
23. Technical Support	13
24. Contact Information	13
25. References	14
26. Symbols	14

1. Introduction

HLA-B27 RQ kit provides a ready-to-use Real-Time polymerase chain reaction (PCR) test designed for detecting HLA-B27 sequence. All required reagents are included in the PCR Mix provided in the kit. This Mix also contains a different series of primers and probe for detecting a housekeeping gene sequence as an Internal Control (IC). IC prevents false negative results due to extraction or presence of PCR inhibitors.

This kit is intended for Research Use Only!

2. Intended Use

HLA- B27 RQ kit is intended for detecting HLA-B27 in human DNA. Detection is achieved using Real-Time PCR and is compatible with Rotor-Gene, StepOne and MIC machines.

3. Background Information

Human leukocyte antigen (HLA-B27) is a class I surface antigen encoded by the B locus in the major histocompatibility complex (MHC) on chromosome 6. HLA-B27 has been associated with ankylosing spondylitis and other inflammatory diseases such as reactive arthritis, uveitis, and psoriasis. Studies show 90% of patients with ankylosing spondylitis have HLA-B27 compared with 5-10% of healthy controls. As a result, tests for HLA-B27 are valuable in the diagnosis and control of these diseases.

This kit detects 98.5% of HLA-B27 subtypes listed in the IMGT/HLA (Release 3.50.0 (2022-10-12)). All 389 subtypes of HLA-B27 detected with this kit except B27:18, B27:23, B27:29, B27:75, B27:157, and B27:189. This kit also detects 6 non-B27 alleles including B38:22, B40:75, B40:322, B44:97, B44:263, B49:26 which results in a specificity of 98%.

4. Test Principle

The target sequence is detected using PCR, where primers specific to the target genome amplify its unique sequence. Real-Time PCR facilitates the detection of the amplified product through fluorescent-labeled probes. Therefore, monitoring fluorescence provides a means for detecting the target without requiring post-amplification analysis. This eliminates the possibility of PCR product contamination.

5. Kit Contents

The kit contains a manual, a flash card, and the following reagents:

Label	Content	Quantity
HLA-B27 RQ Mix	PCR Mix*	480 µl
Pos Ctrl	Positive Control	100 µl
Neg Ctrl	Negative Control	100 µl
Water	PCR Grade Water	200 µl

*1, 2 or 4 tubes for 24, 48 or 96 reaction kits.

6. Packaging models

The kit is available in 24, 48, and 96 reactions of 25 microliters.

7. Storage and Stability

The kit components should be shipped and stored at -20°C and are stable until the expiration date mentioned. Avoid repeated freeze-thaw more than three times to prevent reduced sensitivity.

8. Product Use Limitations

- This kit is intended to be used only by specially instructed and trained personnel.

- The User manual should be strictly followed, and any modification will invalidate the results.
- The kit and its contents should not be used past the expiration date on the package.
- The kit and its contents should not be used if there is any sign of pink or red color on the Warm Mark label.
- This kit is for Research use only and is not validated for IVD (in vitro diagnostics) applications.

9. Additionally Required Materials

To use this kit, you need the following items:

- Real-Time PCR machine and accessory computer
- Tabletop microtube centrifuge
- Vortex Mixer
- Dry Block Heater
- Adjustable pipettors and nuclease free filtered tips
- DNA extraction kit
- Nuclease free 1.5ml microtubes and PCR microtubes
- Disposable powder-free gloves
- Cold block

10. General Precautions

To prevent false results, always pay attention to the following points:

- **Treat all samples as potentially infectious.**
- Within the pre-PCR work area, assign three separate spaces for:
a) Sample storage and extraction, b) Reagent preparation where the PCR mix is aliquoted into tubes, and c) Reaction preparation area for addition of templates to the tubes.
- Always wipe the working surfaces with 70% Ethanol before and after work.

- **Thaw kit components on crushed ice** completely, mix by flickering followed by a quick spin and **store on crushed ice after**.
- Keep PCR Mix at -20°C at all times. Take it out just before use and return it to freezer immediately after.
- Do not place PCR tubes on crushed ice. Use cold blocks instead.

11. Specimen, Storage and Transport

Proper sample to test is peripheral blood. Whole blood should be collected in sterile condition in proper and sterile tubes. We recommend EDTA or Citrate as anticoagulant. Samples should be shipped and stored at +4°C (stable for a 48 hours). For longer terms, sample should be aliquoted and stored at -20°C which is stable for a few months.

12. Interfering Substances

Heparin (more than 10 IU/ml) affects the PCR. Blood collected in heparin containing tubes and samples of heparinized patients must not be used.

Elevated levels of bilirubin (≤ 4.5 mg/dl) and lipids (≤ 1000 mg/dl) and hemolytic samples do not influence the extraction and PCR.

13. Internal Control

To examine DNA extraction quality as well as the presence of PCR inhibitors and to prevent false negative results, primers, and probe for an Internal Control (a housekeeping gene) are included in the PCR Master Mix. Internal control should generate a CT of 20-35. See analysis section for more details.

14. DNA Isolation

DNA isolation can be performed using different kits from various

manufacturers. We recommend using following:

- High Pure PCR Template Preparation Kit (Cat. no. 11796828001, Roche Applied Science, Mannheim, Germany).
- QIAamp DNA Blood Mini Kit (Cat. no. 51104, Qiagen GmbH, Hilden, Germany).

15. PCR Protocol

Thaw the reagents on crushed ice completely, followed by a brief mixing and a quick spin. Place the required number of microtubes on a cold block. Consider one tube for each sample plus one for each control and one for the NTC.

Pipette 20µl of HLA-B27 RQ Mix directly to each tube followed by adding 5ul of controls, or sample DNA.

Cap the tubes and visually inspect to make ensure all are capped securely. Place the tubes in the machine.

Note: Working with StepOne instrument, spin tubes briefly before loading on the block.

Note: If using Rotor-Gene attach the locking ring too.

16. Devices and software

HLA-B27 RQ kit is designed to work with Rotor-Gene, StepOne and MIC.


17. Programming Rotor-Gene

- Before you start the machine, make sure you have attached the locking ring on the rotor!

Open the HLA-B27 template file for Rotor-Gene (provided in the flash card, or accessible by kit QR code); HLA-B27 0.1 is for strip tubes and HLA-B27 0.2 is for 0.2ml tubes. Program starts.

Note: For Gain Optimisation, in the View menu select the Gain Optimisation. Adjust the setting according to the below image.

Make sure to set the Tube Position to number 1 for all channels (note that Tube number 1 should contain HLA-B27 Mix).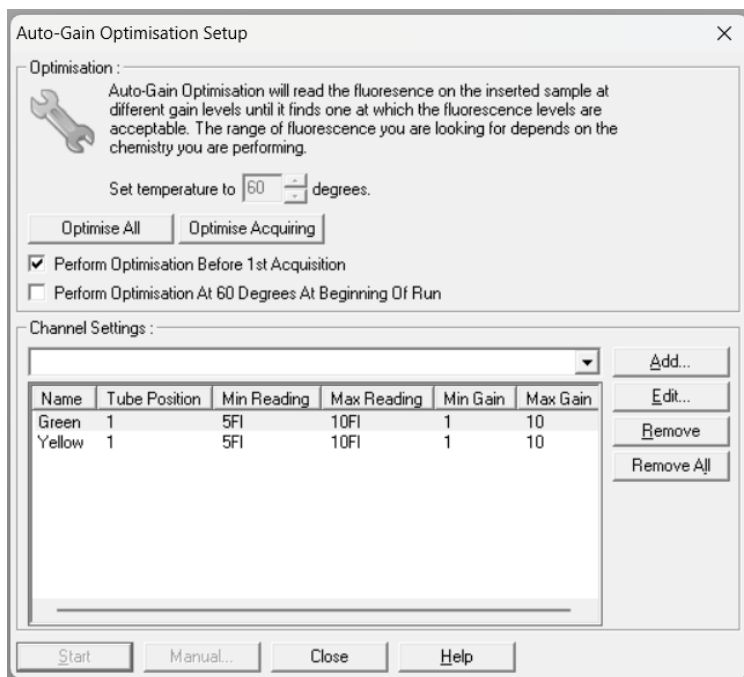

Click on Start button (Green button on the top menu). On the pop-up window click Start again and save the run file. Edit sample names.

18. Programming StepOne

Open the StepOne software (V 2.*). On the Set Up menu, click on Template (provided in the flash card, or accessible by kit QR code). Click on "Plate Setup". Controls and few samples are defined. You may change plate setup using right-click options (copy, paste, clear). You may also add or remove samples on "Define Targets and Samples" menu. When finished, click on "Start Run" and save the experiment to the desired location. Instrument will start shortly.

19. Programming Other Machines

If you apply this kit to other Real-Time PCR machines, program it according to the following table:

Step	Temperature and time	Cycles
1	95°C x 3 min	1
2	95°C x 15 sec	45
	60°C x 60 sec	

Fluorescence should be collected at 60°C for FAM and VIC dyes. PCR Mix contains ROX with the final concentration of 300nM in the reaction.

20. Data Analysis: Rotor-Gene

Before analyzing results, make sure that in the sample menu, Patient samples are defined as "unknown" and Negative control or no template control as "Negative Control" or "NTC" respectively. To analyze data briefly, click on the Analysis menu and then under the Quantitation tab double-click on Cycling A. Green. Manually put threshold at 0.1. To analyze data briefly, click on the analysis menu and then under the Quantitation tab double-click on cycling A. Green. Put threshold at 0.1 Manually. Repeat the above for the Yellow Channel. Figures 1 and 2 represent typical graphs for the Rotor-Gene. To interpret the results, please note that a signal in the **Green** channel is due to **HLA-B27**, and a signal in the **Yellow** channel is due to the **IC**.

Note that a sample is considered Positive only if it has a sigmoid graph and log phase, and only then CT is reliable and can be used. In the absence of a sigmoid graph and log phase, the sample is considered Negative, and CT, if present, is not reliable.

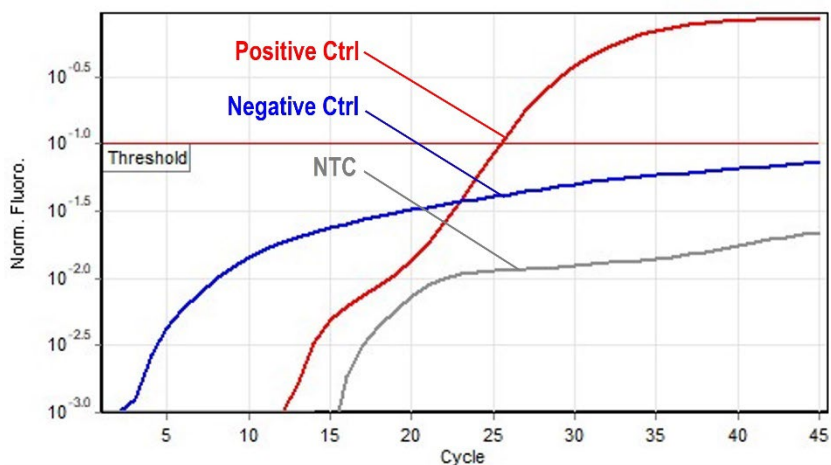
Fig 1. Typical HLA-B27 graph in Green channel for Rotor-Gene

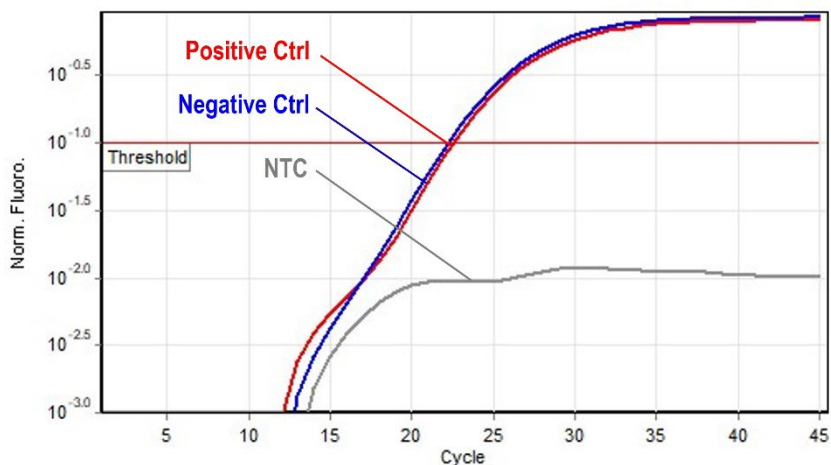
Fig 2. Typical HLA-B27 graph in Yellow channel for Rotor-Gene

Consider the following points when analyzing:

- If a sample is positive in both Green and the Yellow channels with sigmoid graphs and delta-CT between the two channels is less than 5 cycles, then the sample is **Positive** for HLA-B27.
- A sample is **Negative** for HLA-B27 if it is negative in the Green channel while it is positive for Yellow channel with a sigmoid graph and CT of 20-35.
- A sample is **Negative** for HLA-B27 if it is positive in the Green channel with CT of 25-40 while it is positive for Yellow channel with a sigmoid graph and CT of 20-35 and delta-CT between Green and Yellow is more than 5 cycles.
- Results are **Inconclusive** and the test should be repeated if a sample is negative in the Yellow channels.

The interpretation of results is summarized in the following Table.

Δ CT: Green – Yellow	Green/FAM	Yellow/VIC	Result
Δ CT < 5 in RG Δ CT < 5 in StepOne	+ CT: 20-40	+ CT: 20-35	Positive
Δ CT > 5 in RG Δ CT > 5 in StepOne	+ CT: 25-40	+ CT: 20-35	Negative
–	–	+ CT: 20-35	Negative
–	– / +	–	Inconclusive

21. Data Analysis: StepOne

Analyze data according to StepOne manual. Briefly, click on Analyze and set the threshold for the **FAM** and for the **VIC** at 0.1. Please note that signal in the **FAM** channel is due to **HLA-B27**, and a signal in the **VIC** channel is due to the **IC**.

Note that a sample is considered Positive only if it has a sigmoid graph and log phase, and only then CT is reliable and can be used. In the absence of a sigmoid graph and log phase, the sample is considered Negative, and CT, if present, is not reliable.

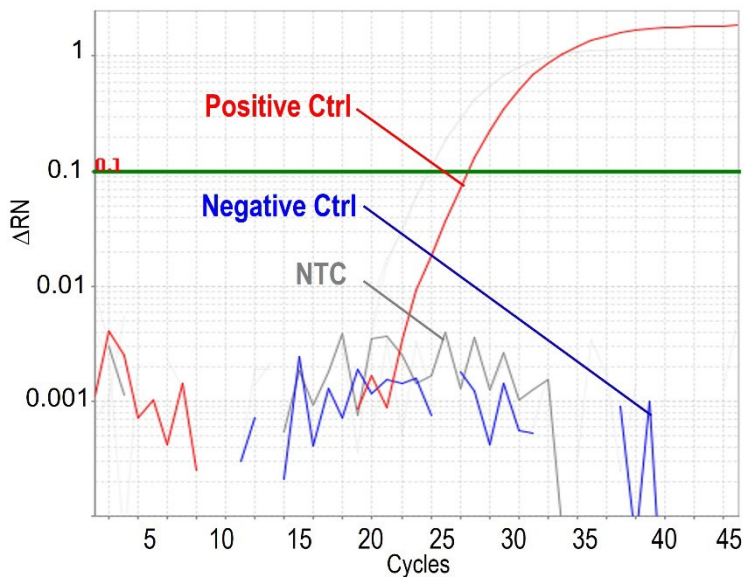
Fig 3. Typical HLA-B27 graph in FAM channel for StepOne

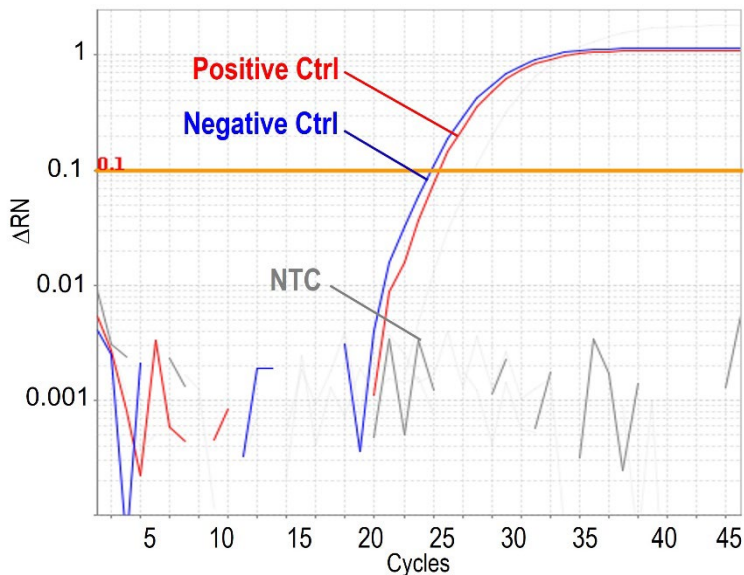
Fig 4. Typical HLA-B27 graph in VIC channel for StepOne

Consider the following points when analyzing:

- If a sample is positive in both the FAM and the VIC channels with sigmoid graphs and delta-CT between the channels is less than 5 cycles, then the sample is **Positive** for HLA-B27.
- A sample is **Negative** for HLA-B27 if it is negative in the FAM channel while it is positive for VIC channel with a sigmoid graph and CT of 20-35.
- A sample is **Negative** for HLA-B27 if it is positive in the FAM channel with CT of 25-40 while it is positive for the VIC channel with a sigmoid graph and CT of 20-35 and delta-CT between FAM and VIC is more than 5 cycles.
- Results are **Inconclusive** and the test should be repeated if a sample is negative in VIC channel.

Interpretation of results is summarized in the page 10, Table.

22. Disposal Method

The contents of the kit do not require any special treatment before disposal and can be directly discarded. Infectious specimens should be maintained in 5% Sodium Hypochlorite overnight and then discarded.

23. Technical Support

For technical support, contact us via
Phone at +98-9936223241
email: info@novingene.com

24. Contact Information

NovinGene ParsVira

Address: No. 5, Iranshahr St, Tehran, Iran 1581633336.

Tel: +98 21-88837393

+98 990-1813124

Email: info@novingene.com

Website: www.novingene.com

25. References

- Fernando, M.M., Stevens, C.R., Walsh, E.C., De Jager, P.L., Goyette, P., Plenge, R.M., Vyse, T.J. and Rioux, J.D., 2008. Defining the role of the MHC in autoimmunity: a review and pooled analysis. PLoS genetics, 4(4), p.e1000024.
- Jung, J.H., Bang, C.H., Seok, H., Choi, S.J. and Song, G.G., 2019. Clinical findings of ankylosing spondylitis with and without Human Leukocyte Antigen (HLA)-B27 and HLA-B51. Ann Acad Med Singapore, 48(10), pp.321-9.
- Mackay IM. Real-time PCR in the microbiology laboratory. Clin. Microbiol. Infect. 2004; 10 (3): 190 – 212.
- Xi, Y., 2014. HLA and Associated Important Diseases. BoD – Books on Demand.

26. Symbols

RUO Research use only	 Manufacturer	 Consult instructions for use
LOT Lot number	 Content sufficient for <n> tests	 Use-by date
REF Catalogue number	SN Serial number	 Temperature limit

For more information and resources please visit our website; www.novingene.com