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ایمیل: [info@novingene.com](mailto:info@novingene.com)

وبسایت: [www.novingene.ir](http://www.novingene.ir)

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- Alspaugh, J.A., 2015. Virulence mechanisms and *Cryptococcus neoformans* pathogenesis. *Fungal Genetics and Biology*, 78, pp.55-58.
- Bateman, M., Oladele, R. and Kolls, J.K., 2020. Diagnosing *Pneumocystis jirovecii* pneumonia: a review of current methods and novel approaches. *Medical Mycology*, 58(8), pp.1015-1028.
- Fosses Vuong, M., Hollingshead, C.M. and Waymack, J.R., 2023. *Aspergillosis*.
- Gagneux, S., 2018. Ecology and evolution of *Mycobacterium tuberculosis*. *Nature Reviews Microbiology*, 16(4), pp.202-213.
- Mackay IM., 2004. Real-time PCR in the microbiology laboratory. *Clin. Microbiol. Infect.* 10 (3): 190 – 212.
- Narciso, A.R., Dookie, R., Nannapaneni, P., Normark, S. and Henriques-Normark, B., 2025. *Streptococcus pneumoniae* epidemiology, pathogenesis and control. *Nature Reviews Microbiology*, 23(4), pp.256-271.
- Prasad, R., 2017. *Candida albicans*. Springer International Publishing.

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# Bacto-Fungal Respiratory Panel Kit Manual

Spring 2026, Version 1.0

For Real-Time PCR Detection of *Aspergillus* spp., *Candida albicans*,  
*Cryptococcus neoformans*, *Mycobacterium tuberculosis* complex,  
*Pneumocystis jirovecii*, *Streptococcus pneumoniae* DNA  
For Research Use Only

 24 (Cat# RespIIIIRQ24)

 48 (Cat# RespIIIIRQ48)

 96 (Cat# RespIIIIRQ96)


 NG-WI-ASL-74-100

RUO

**NovinGene ParsVira**

No. 20, 4<sup>th</sup> St, Gisha, Tehran, Iran 1446843434.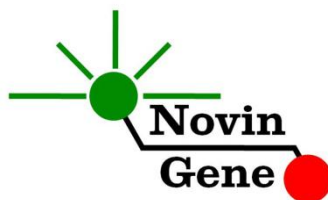

# Table of Contents

1. Introduction .....	3
2. Intended Use .....	3
3. Background Information .....	3
4. Test Principle .....	5
5. Kit Contents .....	5
6. Packaging models .....	6
7. Storage and Stability .....	6
8. Product Use Limitations .....	6
9. Additionally Required Materials .....	6
10. General Precautions .....	7
11. Specimen, Storage and Transport .....	7
12. Internal Control (IC) .....	7
13. DNA Isolation .....	8
14. PCR Protocol .....	8
15. Devices and software .....	9
16. Programming of the Rotor-Gene .....	9
17. Programming Other Machines .....	10
18. Data Analysis: Rotor-Gene .....	11
19. Sensitivity .....	18
20. Disposal Method .....	18

21. Technical Support.....	19
22. Contact Information.....	19
23. References .....	19
24. Symbols.....	20

## 1. Introduction

Bacto-Fungal RQ kit provides a ready-to-use Real-Time PCR test designed for detecting DNA of respiratory pathogens including bacterial pathogens; *Streptococcus pneumoniae* and *Mycobacterium tuberculosis* complex and Fungal pathogens; *Aspergillus* spp., *Candida albicans*, *Cryptococcus neoformans* and *Pneumocystis Jirovecii*. All required reagents are included in the PCR Mixes provided in the kit. This kit also contains a different series of primers and probe for detecting a synthetic DNA sequence. The kit supplies this synthetic DNA sequence as an Internal Control (IC). The IC can be used either during DNA extraction or in the PCR reaction to prevent false negative results due to failure in the above steps.

This kit is intended for Research Use Only!

## 2. Intended Use

The Bacto-Fungal RQ Kit is a multiplex real-time PCR assay designed for the simultaneous detection of six clinically significant bacterial and fungal respiratory pathogens in two separate reactions. These pathogens include, *Mycobacterium tuberculosis* complex, *Streptococcus pneumoniae*, *Aspergillus* spp., *Candida albicans*, *Cryptococcus neoformans* and *Pneumocystis Jirovecii*. This kit is compatible with Rotor-Gene and MIC machines.

## 3. Background Information

*Mycobacterium tuberculosis* is a slow-growing, acid-fast bacillus and the etiologic agent of tuberculosis. Transmission occurs via airborne droplet nuclei, leading primarily to pulmonary infection characterized by granulomatous inflammation and caseating necrosis. It is capable of latent infection with potential for reactivation, particularly under conditions of immunosuppression.

*Streptococcus pneumoniae* is a Gram-positive, encapsulated diplococcus and a major human respiratory pathogen. It is transmitted via respiratory droplets and commonly colonizes the nasopharynx before causing invasive or non-invasive disease. Clinically, it is a leading cause of community-acquired pneumonia, meningitis, otitis media, and bacteremia, particularly in vulnerable populations such as children and the elderly.

*Aspergillus* species are filamentous, saprophytic fungi widely distributed in environmental sources such as soil, air, and decomposing organic matter. They are clinically significant opportunistic pathogens associated with a spectrum of pulmonary diseases, particularly in immunocompromised patients.

*Candida albicans* is an opportunistic dimorphic yeast that is part of the normal human microbiota of mucosal surfaces including the oropharyngeal and gastrointestinal tracts. Although commonly considered a colonizer in respiratory samples, it may be associated with invasive candidiasis in severely immunocompromised patients or those in intensive care settings. Its clinical relevance in respiratory specimens should be interpreted in conjunction with clinical presentation and risk factors.

*Cryptococcus neoformans* is an encapsulated basidiomycetous yeast found in environmental reservoirs such as soil contaminated with avian excreta. It is a clinically important opportunistic pathogen capable of causing pulmonary cryptococcosis following inhalation of infectious propagules. In immunocompromised individuals, causing cryptococcal meningitis.

*Pneumocystis jirovecii* is an atypical fungal organism lacking ergosterol in its cell membrane and it is classified as an obligate extracellular pathogen of the lung alveoli. *P. jirovecii* is associated with *Pneumocystis pneumonia* (PCP), predominantly in immunocompromised hosts such as HIV-infected patients or individuals receiving immunosuppressive therapy.

#### 4. Test Principle

The pathogens are detected using PCR, where primers specific to the target genome amplify its unique sequence. Real-Time PCR facilitates the detection of the amplified product through fluorescent-labeled probes. Monitoring fluorescence provides a means for detecting the target without requiring post-amplification analysis. This eliminates the possibility of PCR product contamination.

#### 5. Kit Contents

The kit contains a manual and the following reagents:

Label	Content	Quantity
Bacto A Mix	PCR mix* for Mycobacterium tuberculosis, Streptococcus pneumoniae, Internal Control	360 µl
Fungal A Mix	PCR mix* for Aspergillus pp., Candida albicans, Cryptococcus neoformans, Pneumocystis Jirovecii	360 µl
Positive Ctrl	Positive Control	250 µl
Internal Ctrl	Internal Control*	250 µl
Water	PCR Grade Water	200 µl

\* 1, 2 and 4 tubes for 24, 48 and 96 reaction kits respectively.

## 6. Packaging models

The kit is available in 24, 48, and 96 reactions of 25 microliters.

## 7. Storage and Stability

The kit components should be shipped and stored at  $-20^{\circ}\text{C}$  and are stable until the expiration date mentioned. Avoid repeated freeze-thaw more than three times to prevent reduced sensitivity.

## 8. Product Use Limitations

- This kit is intended to be used only by specially instructed and trained personnel.
- The user manual should be strictly followed, and any modification will invalidate the results.
- The kit and its contents should not be used past the expiration date on the package.
- The kit and its contents should not be used if there is any sign of pink or red color on the Warm Mark label.
- This kit is for Research use only and is not validated for IVD (in vitro diagnostics) applications.

## 9. Additionally Required Materials

To use this kit, you need the following items:

- Real-Time PCR machine and the and the accessory computer
- Tabletop microtube centrifuge
- Vortex Mixer
- Adjustable pipettors and nuclease free filtered tips
- DNA extraction kit and required equipments/items
- PCR microtubes
- Disposable powder-free gloves

- Cold block

## 10. General Precautions

In order to prevent false results, always pay attention to the following points:

- **Treat all samples as potentially infectious. respiratory samples may contain airborne pathogens and should be processed using appropriate containment procedures and personal protective equipment.**
- Within the pre-PCR work area, assign three separate spaces for: a) Sample storage and extraction, b) Reagent preparation where the master-mix is aliquoted into tubes, and c) Reaction preparation area for addition of extracted DNA to the PCR tubes.
- Always wipe the working surfaces with 70% Ethanol before and after work.
- **Place kit components on ice till completely thawed, mix by flickering followed by a quick spin and store on crushed ice after.**
- **Keep the PCR Mix tube at -20°C at all times. Take it out just before use and return it to freezer immediately after.**
- Do not place 0.2ml PCR tubes on crushed ice. Use cold block instead.

## 11. Specimen, Storage and Transport

Samples should be collected from lower respiratory tract and include bronchoalveolar lavage (BAL) and Sputum. Samples could be stored at 2-8°C for a few hours and at -20°C or lower for longer periods.

## 12. Internal Control (IC)

To assess the possibility of DNA extraction failure and PCR inhibition and to prevent false negative results, this kit contains an

internal control (IC). IC should be used during the extraction process to monitor both DNA extraction and PCR reaction. IC should be added to the mixture of lysis buffer and patient sample during extraction. The required volume of IC is 10% of the elution buffer. For instance, if the extracted DNA is eluted with 100ul, then 10ul of IC should be added to the mixture of lysis buffer and patient sample. **Please note that IC should not be added directly to the patient sample (i.e. before the addition of lysis buffer) as it loses its efficiency.**

In a successful DNA extraction and PCR test, IC should generate a CT of 28-32 in Yellow channel with Mix Bacto A.

### 13. DNA Isolation

DNA isolation can be performed using different kits from various manufacturers. We recommend the following:

- High Pure PCR Template Preparation Kit (Cat# 11796828001, Roche Applied Science, Mannheim, Germany)

### 14. PCR Protocol

This kit includes two mixes, **Bacto A Mix** which detects *Streptococcus pneumoniae* and *Mycobacterium tuberculosis* complex and **Fungal A Mix** which detects *Aspergillus* spp., *Candida albicans*, *Cryptococcus neoformans*, and *Pneumocystis Jirovecii*.

Thaw the reagents on ice completely followed by a brief mixing and a quick spin.

Each sample should be examined in two reactions in parallel. Place two series of tubes on cold block. In each series, consider one tube for each sample plus one for positive control and one for water.

**Aliquot 15ul of Bacto A Mix to the first series of PCR tubes**

**Aliquot 15µl of Fungal A Mix to each tube of the second series followed by adding 10µl of extracted DNA, Positive Control or water.**

Cap the tubes and visually inspect to ensure all are capped securely. Place the tubes in the machine.

*Note: If using Rotor-Gene attach the locking ring.*

## **15. Devices and software**

Bacto-Fungal RQ kit is designed for use with Rotor-Gene and MIC.


## **16. Programming of the Rotor-Gene**

*Before starting the machine, ensure you have attached the locking ring on the rotor!*

Open the Bacto-Fungal template file for Rotor-Gene (provided in the Flash card, or accessible by Kit QR code) double-click on "Bacto-Fungal 0.1" or " Bacto-Fungal 0.2" depending on the tubes used.

Note: in the View menu, select the Gain Optimisation. Under the Auto-Gain Optimisation Setup, select the Optimise Acquiring option. Adjust the setting according to the following image.

make sure the selected Tube Position contains Fungal A Mix. Enable the option "Perform Optimisation Before 1st Acquisition" and close the window.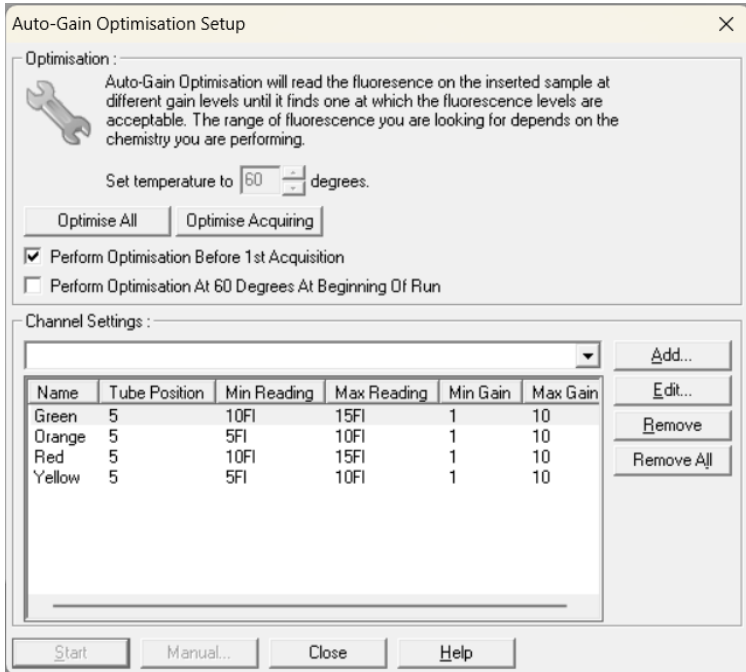

Gain Optimisation Setting

Click on the Start button (Green button on the top menu). On the pop-up window click Start again and save the program.

## 17. Programming Other Machines

If you apply this kit to other Real-Time PCR machines, program it according to the following table:

Step	Temperature and time	Cycles
1	<b>95°C x 3 min</b>	1
2	<b>95°C x 15 sec</b>	45
	<b>60°C x 75 sec*</b>	

Fluorescence should be collected at 60°C for FAM, VIC, ROX and Cy5 dyes. *In this experiment, ROX should not be selected as the reference dye.*

\* 60 seconds for Rotor-Gene, 75 seconds for other machines.

## 18. Data Analysis: Rotor-Gene

The results are interpreted based on the reaction of the samples with the two mixes according to the following table.

Mix/Channel	Green/FAM	Yellow/HEX	Orange/ROX	Red/Cy5
Bacto A Mix	Mycobacterium tuberculosis	Internal Ctrl	-	Streptococcus pneumoniae
Fungal A Mix	Cryptococcus neoformans	Pneumocystis Jirovecii	Candida albicans	Aspergillus spp.

### A) Analysis of the sample results with Bacto A Mix:


To analyze data of **Bacto A Mix** briefly, click on the analysis menu and then under Quantitation tab double click on cycling A. Green. Manually put threshold, at 0.1. Repeat the above for Red and Yellow Channels.

Figures 1, 2 and 3 represent typical graphs with Bacto A Mix for Rotor-Gene.


To interpret the results, please note that:

with the **Bacto A Mix**, a signal detected in the **Green** channel indicates **Streptococcus pneumoniae**, a signal in the **Red** channel represents **Streptococcus pneumoniae** and a signal in the **Yellow** channel is due to **Internal Control**.


**Note that a sample is considered Positive only if it has a sigmoid graph and log phase, and only then CT is reliable and can be used. In the absence of a sigmoid graph and log phase, the sample is considered Negative, and CT, if present, is not reliable.**

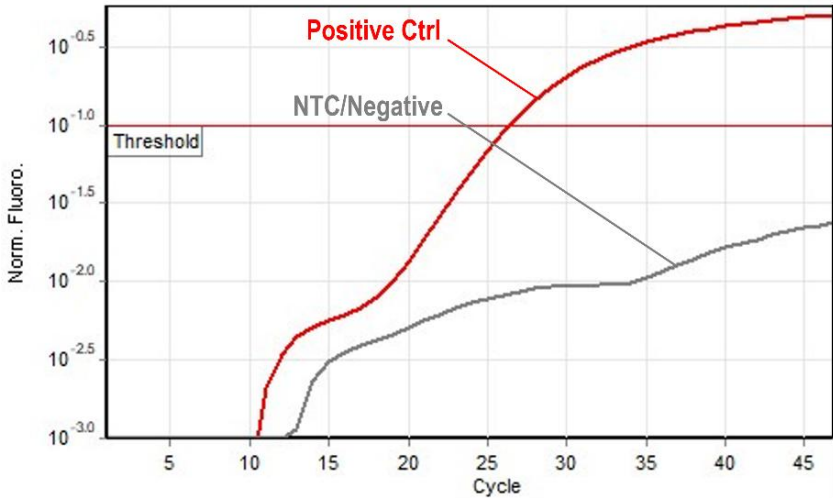**Fig 1.** Typical Controls graph with Bacto A Mix in Green channel for Rotor-Gene (*Mycobacterium tuberculosis* complex)

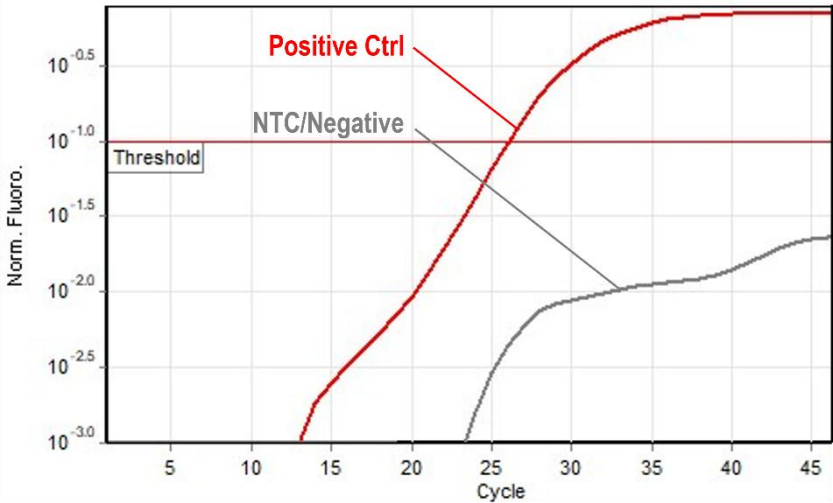**Fig 2.** Typical Controls graph with Bacto A Mix in Red channel for Rotor-Gene (*Streptococcus pneumoniae*)

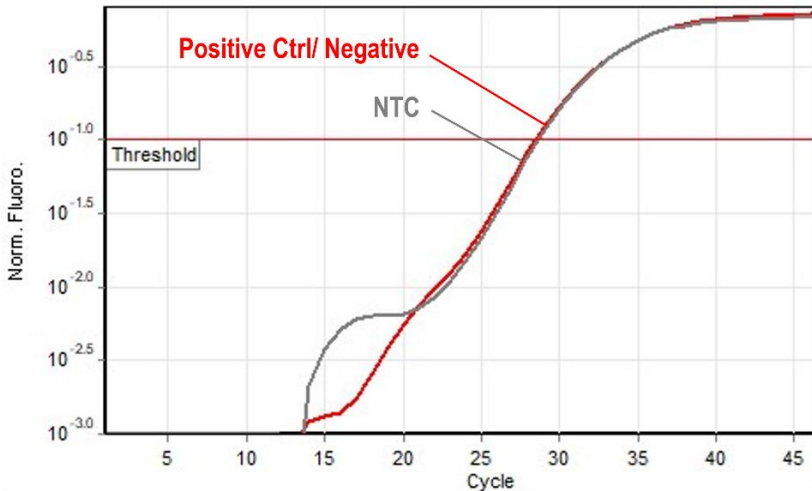**Fig 3.** Typical Controls graph with Bacto Mix in Yellow channel for Rotor-Gene (Internal Control)

Consider the following points when analyzing results with [Bacto A Mix](#):

- A sample is **Positive** for Mycobacterium tuberculosis complex if it is positive in the **Green** channel with a sigmoid graph and CT of less than 40.
- A sample is **Positive** for Streptococcus pneumoniae if it is positive in the **Red** channel with a sigmoid graph and CT of less than 40.
- A sample is **Negative** for Mycobacterium tuberculosis complex and Streptococcus pneumoniae if it is negative in the **Green** and the **Red** channels while it is positive in the **Yellow** channel with a sigmoid graph and CT of 28-32.
- Results are **Inconclusive** and the test should be repeated if a sample is negative in all three channels, **Green**, **Red** and **Yellow** channel with Bacto A Mix.

The interpretation of results with **Bacto A Mix** is summarized in the following Table.


Bacto A Mix			
Green/ FAM	Yellow/ VIC	Red/ Cy5	Result
+	-/+	-	Pos for Mycobacterium tuberculosis
-	-/+	+	Pos for Streptococcus pneumoniae
-	+	-	Negative
-	-	-	Inconclusive

**B) Analysis of the sample results with Fungal A Mix:**


To analyze data of **Fungal A Mix** briefly, click on the analysis menu and then under Quantitation tab double click on cycling A. Green. Manually put threshold, at 0.1 for Green, Orange, Red and Yellow channels.

Figures 4, 5, 6 and 7 represent typical graphs with Fungal A Mix for Rotor-Gene.


A signal with **Fungal A Mix** in the **Green** channel indicates **Cryptococcus neoformans**, in the **Orange** channel indicates **Candida albicans** and in the **Red** channel indicates **Aspergillus spp.** and, in the **Yellow** channel indicates **Pneumocystis Jirovecii**.

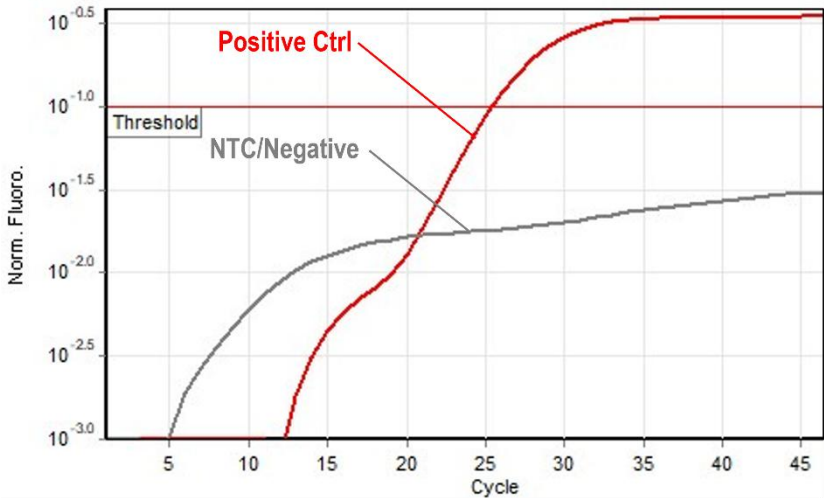
**Fig 4.** Typical Controls graph with Fungal A Mix in Green channel for Rotor-Gene (*Cryptococcus neoformans*)

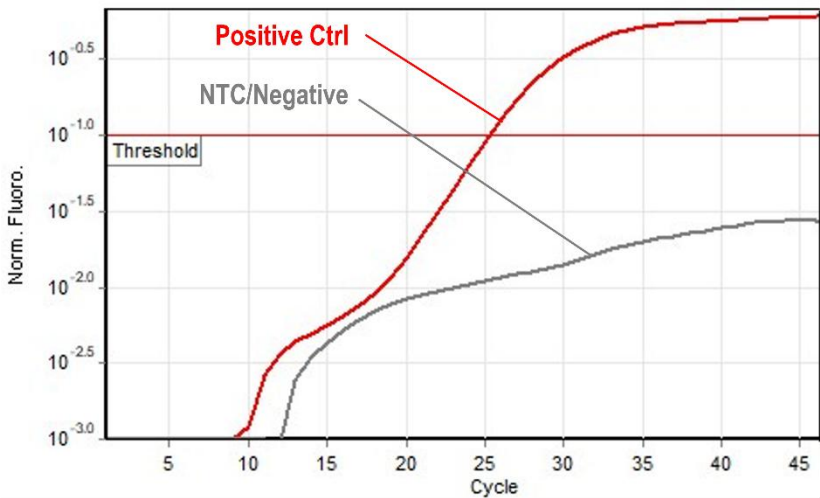**Fig 5.** Typical Controls graph with Fungal A in Yellow channel for Rotor-Gene (*Pneumocystis Jirovecii*)

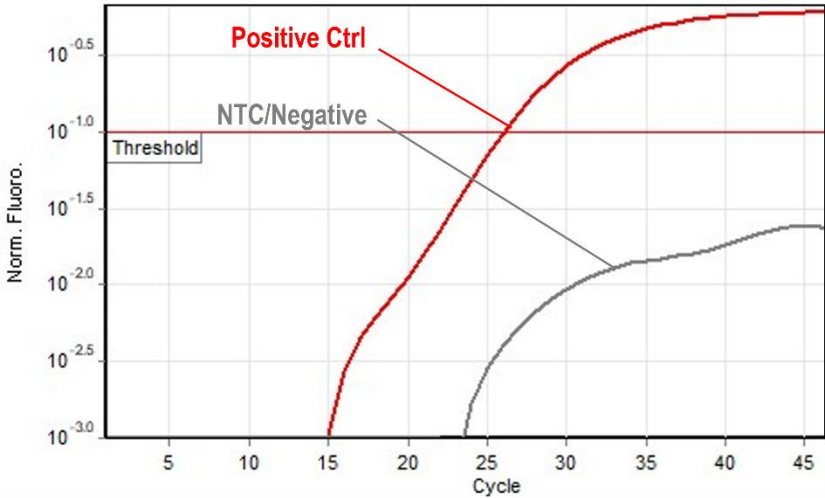**Fig 6.** Typical Controls graph with Fungal A Mix in Red channel for Rotor-Gene (*Aspergillus Spp*)

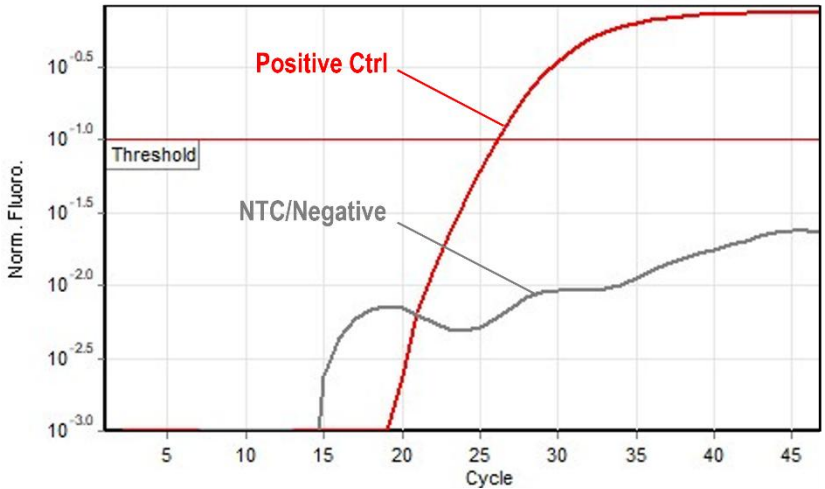**Fig 7.** Typical Controls graph with Fungal A Mix in Orange channel for Rotor-Gene (*Candida albicans*)

**Note that a sample is considered Positive only if it has a sigmoid graph and log phase, and only then CT is reliable and can be used. In the absence of a sigmoid graph and log phase, the sample is considered Negative, and CT, if present, is not reliable**

Consider the following points when analyzing results with **Fungal A Mix:**

- A sample is **Positive** for Cryptococcus neoformans if it is positive in the **Green** channel with sigmoid graphs and CT of less than 40.
- A sample is **Positive** for Candida albicans if it is positive with Fungal A Mix in the **Orange** channel with sigmoid graphs and CT of less than 40.
- A sample is **Positive** for Aspergillus spp. if it is positive in the **Red** channel with sigmoid graphs and CT of less than 40.
- A sample is **Positive** for Pneumocystis Jirovecii if it is positive in the **Yellow** channel with sigmoid graphs and CT of less than 40.
- A sample is **Negative** for Cryptococcus neoformans, Candida albicans, Aspergillus spp. and Pneumocystis Jirovecii if it is negative in the **Green**, the **Orange**, the **Red** and the **Yellow** channels with Fungal A Mix while IC is positive with Bacto A Mix in the **Yellow** channel with a sigmoid graph and CT of 28-32.
- Results are **Inconclusive** and the test should be repeated if a sample is negative with in all four channels with Fungal A Mix and Bacto A Mix.

The interpretation of results for Fungal A Mix is summarized in the following Table.

Fungal A Mix				
Green/ FAM	Yellow/ VIC	Red/ Cy5	Orange/ ROX	Result
+	-	-	-	Positive for <b>Cryptococcus neoformans</b>
-	+	-	-	Positive for <b>Pneumocystis Jirovecii</b>
-	-	+	-	Positive for <b>Aspergillus spp.</b>
-	-	-	+	Positive for <b>Candida albicans</b>
-	-	-	-	Negative if sample IC is <u>Positive</u>
-	-	-	-	Inconclusive (If sample is also Negative in all channels with Bacto A Mix)

### 19. Sensitivity

The analytical detection limit of the kit was assessed with the cloned targets genomes (partial) and showed a limit of detection equal to 1 copies/μl for Mycobacterium tuberculosis, 38 copies/μl for Streptococcus pneumoniae, 18 copies/μl for Cryptococcus neoformans, 87 copies/μl for Aspergillus spp., 21 copies/μl for Pneumocystis Jirovecii and 24 copies/μl for Candida albicans.

### 20. Disposal Method

The contents of the kit do not require any special treatment before disposal and can be directly discarded. Infectious specimens should be maintained in 5% Sodium Hypochlorite overnight and

then discarded.

## 21. Technical Support

For technical support, contact us via

Phone: +98 993-6223241

Email: [info@novingene.com](mailto:info@novingene.com)

## 22. Contact Information

### NovinGene ParsVira

Address: No. 20, 4<sup>th</sup> St, Gisha St, Tehran, Iran 1446843434.

Tel: +98 21-88837393

+98 990 -1813124

Email: [info@novingene.com](mailto:info@novingene.com)

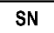
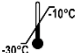
Website: [www.novingene.com](http://www.novingene.com)

## 23. References

- Alspaugh, J.A., 2015. Virulence mechanisms and *Cryptococcus neoformans* pathogenesis. *Fungal Genetics and Biology*, 78, pp.55-58.
- Bateman, M., Oladele, R. and Kolls, J.K., 2020. Diagnosing *Pneumocystis jirovecii* pneumonia: a review of current methods and novel approaches. *Medical Mycology*, 58(8), pp.1015-1028.
- Fosses Vuong, M., Hollingshead, C.M. and Waymack, J.R., 2023. Aspergillosis. "
- Gagneux, S., 2018. Ecology and evolution of *Mycobacterium tuberculosis*. *Nature Reviews Microbiology*, 16(4), pp.202-213.
- Mackay IM., 2004. Real-time PCR in the microbiology laboratory. *Clin. Microbiol. Infect.* 10 (3): 190 – 212.

- Narciso, A.R., Dookie, R., Nannapaneni, P., Normark, S. and Henriques-Normark, B., 2025. Streptococcus pneumoniae epidemiology, pathogenesis and control. Nature Reviews Microbiology, 23(4), pp.256-271.
- Prasad, R., 2017. Candida albicans. Springer International Publishing.

## 24. Symbols

 <b>RUO</b> Research use only	 <b>Manufacturer</b>	 <b>Consult instructions for use</b>
 <b>LOT</b> Lot number	 <b>Content sufficient for &lt;n&gt; tests</b>	 <b>Use-by date</b>
 <b>REF</b> Catalogue number	 <b>SN</b> Serial number	 <b>Temperature limit</b>

**For more information and resources please visit our website; [www.novingene.com](http://www.novingene.com)**